

PHYC 500: Introduction to LabView

M.P. Hasselbeck, University of New Mexico

Exercise 9 (v 1.1) Spectral content of waveforms

This exercise provides additional experience with the Waveform palette, along with more cluster and array operations. A VI will be setup to produce a modulated sine wave followed by Fourier analysis.


Open a blank VI and go to the Block Diagram. Right-click and select Waveform: Analog Wfm: Generation: Sine Wfm. Alternatively, locate it with Quick Drop. Place two of these icons on the Block Diagram. One will generate the fundamental or carrier wave at frequency f_0 and the second will be the modulation signal at frequency f_1 . Note that by definition $f_0 > f_1$. Create a Front Panel numerical control on the **frequency** terminal of the first sine-wave generator and call it “Fundamental frequency (f_0)”. Do the same on the second and label it “Modulation frequency (f_1)”. Create a control on one generator for the terminal **sampling info**; this will produce a brown cluster icon on the Block Diagram. Wire this to the second Sine Wfm VI so that both are setup the same way. This control cluster sets the waveform sampling rate (Fs) and number of points (#s) in the signal. The defaults are 1 kHz and 1000 points, respectively.

Next multiply the **signal out** terminals together using the multiplication function on the Numeric Palette. Create an indicator on the Multiply output terminal; it will be a waveform cluster. Double-click on the icon to locate the indicator on the Front Panel. Right-click on the edge of the cluster and Replace: Graph: Waveform Graph. It is important to select the entire cluster and not one of its components (eg. **t0**, **dt**, or **Y**). This graph will be the modulated output waveform.


In addition to the output waveform, both input signals should be monitored. Select Waveform: Get Waveform Components and place on the Block Diagram. Wire the “signal out” from the Sine Wave generator at f_0 to the input. Click and drag the components icon down to show two output terminals (**Y** and **attributes**). Alternatively, right-click and “Add Element”. Right-click on the lower terminal, Select Item: dt. This corresponds to the time separation between points. Wire a second “Get Wfm Components” VI to the **signal out** terminal of the second Sine Wave generator that produces frequency f_1 . Only the Y data is needed here.

The two waveforms will be plotted on the same Waveform Graph. In this exercise, the needed information is bundled into a cluster. This cluster contains the starting x -value, dt, and a 2-D array with data from both waveforms. Select Cluster, Class, & Variant:

Bundle and place it on the Block Diagram. Right-click on the top input terminal and create a floating point (DBL) constant of 0. Wire the second input terminal to the dt value from the first waveform. The lower terminal will connect to a 2-D array with both sets of y-data. Select Array: Build Array and re-size it to have two inputs. Connect the 1-D arrays (indicated by Y) to the two inputs of Build Array and wire the 2-D output (double orange wire) to the bottom terminal of the Bundle function. Consult the example Block Diagram below.

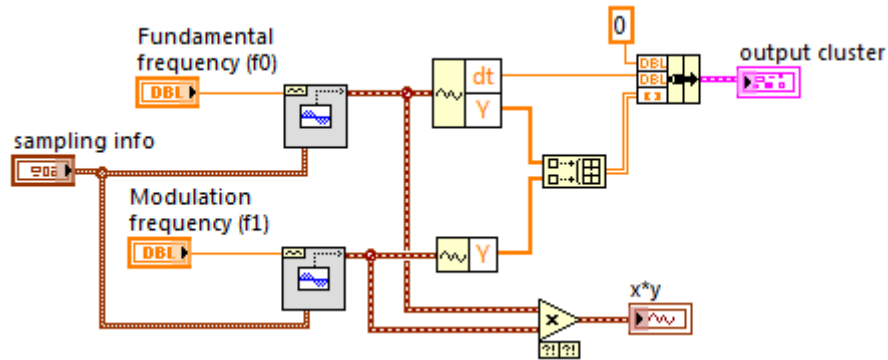

Create an indicator at the output of the Bundle operation; double-click on the icon to locate it on the Front Panel. Right-click on the edge of the cluster and Replace: Graph: Waveform Graph. As before, be careful to click on the cluster boundary and not the internal components. You should now have a working VI. Select a modulation frequency smaller than the fundamental frequency ($f_1 < f_0$) and verify operation. An example Front Panel is shown below.

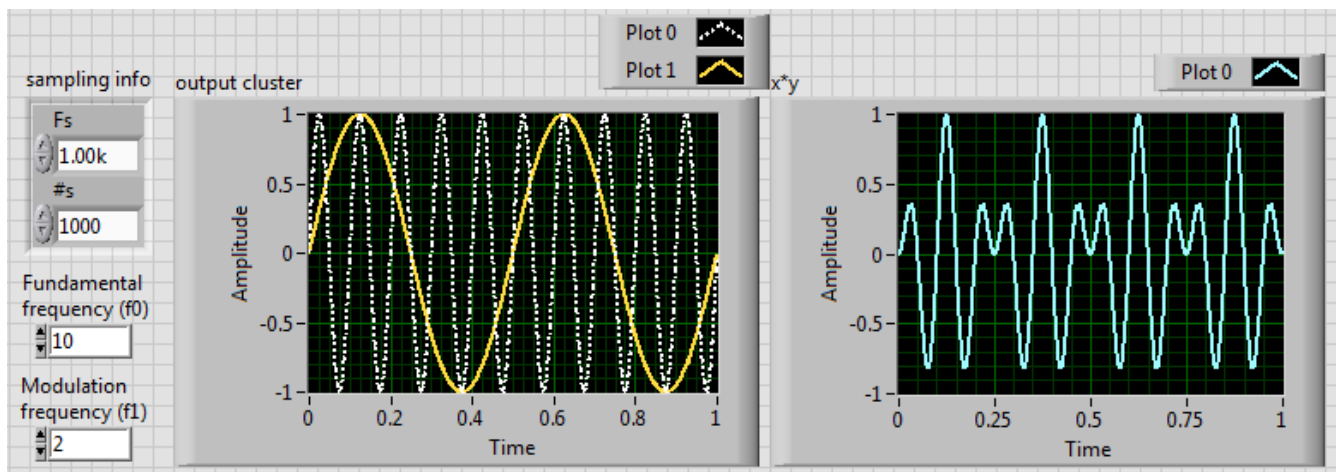


The waveforms on the graphs can be easily customized. Drag down the Plot Legend on the multi-plot graph to reveal the second channel and right-click.

Fourier Analysis

The Fourier power spectrum of the modulated signal is given by the squared amplitude of its Fourier transform $|F(v)|^2$. Select Waveform: Analog Wfm: Measurements: FFT Power Spectrum and PSD.vi and place it on the Block Diagram. Connect a wire from the modulated waveform to the **time signal** input terminal. Create an indicator on the **Power Spectrum/PSD** output terminal. Double-click on the icon to find it on the Front Panel and replace the cluster with a Waveform Graph.

Run the VI and examine the FFT graph. Right-click on the Plot Legend icon and select the line and symbol plot. You should see two peaks at frequencies $f = f_0 \pm f_1$. The sampling frequency F_s will set the time step dt of the temporal waveform, which defines the maximum frequency in the Fourier spectrum. Choice of these parameters may make it very difficult to resolve peaks in the power spectrum graph. One option is to zoom-in on the graph. Right-click on the graph and select Visible Items: Graph Palette. Select the middle tool in the top row to re-size the x -axis display. The entire display can be restored by clicking on the first tool in the second row.

Another option is to edit the displayed data range. Right-click on the graph, select Properties: Scales. Disable Autoscale and set an appropriate Maximum for the x -axis. This operation can also be performed directly on the right-most scale marker on the x -axis. Simply select it and change its value. An FFT power spectrum for $f_0 = 10$ Hz, $f_1 = 2$ Hz, and $f_{\max} = 20$ Hz is shown below.

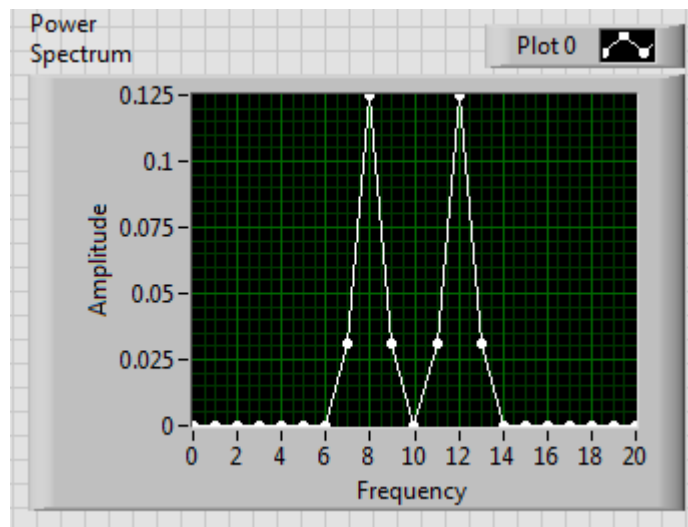


The maximum value of the x -scale can also be set programatically inside the VI. For example, the upper-limit of the frequency axis can be set to twice the value of f_0 or f_1 , depending on which one is larger. Go to the Block Diagram and select Comparison: Max

and Min. Wire the f_0 and f_1 controls to the two inputs labeled x and y . Multiply the $\max(x,y)$ output by 2. Right-click on the Power Spectrum graph icon and create a Property Node: X Scale: Range: Maximum. Right-click on the icon and select “Change to Write”. Wire a connection from its input terminal to the output of the multiplication operation.


When the VI is run, you will notice that the Property Node fails to limit the maximum value of the frequency axis. This is caused by dataflow execution in the Block Diagram. The code executes as soon as data becomes available; re-arranging the icons will not change the timing. The maximum value on the x -axis is set by the Property Node, but that is overwritten by the FFT VI.

One solution is to force the order of execution with a Flat Sequence structure. Locate this on the Structures palette and place it in the Block Diagram. Drop the Power Spectrum icon inside the structure, right-click on its boundary, and “Add Frame After”. Place the Property Node in the second frame. This rigidly defines the order of execution. The complete Block Diagram will look similar to this:

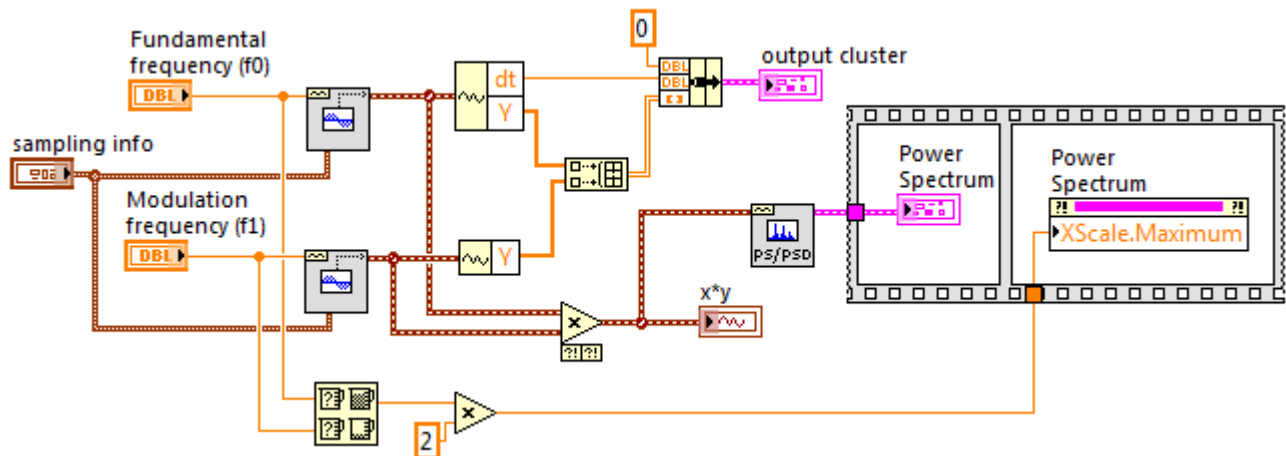

The Flat Sequence first draws the FFT graph, which is immediately followed by an x -axis rescaling.

A simpler and more efficient approach to dataflow programming can sometimes be implemented with error handling terminals. Error information can be forced to propagate through a sequence of component VIs by wiring the error cluster serially in the Block Diagram. This is illustrated with modifications to the above VI as shown below:

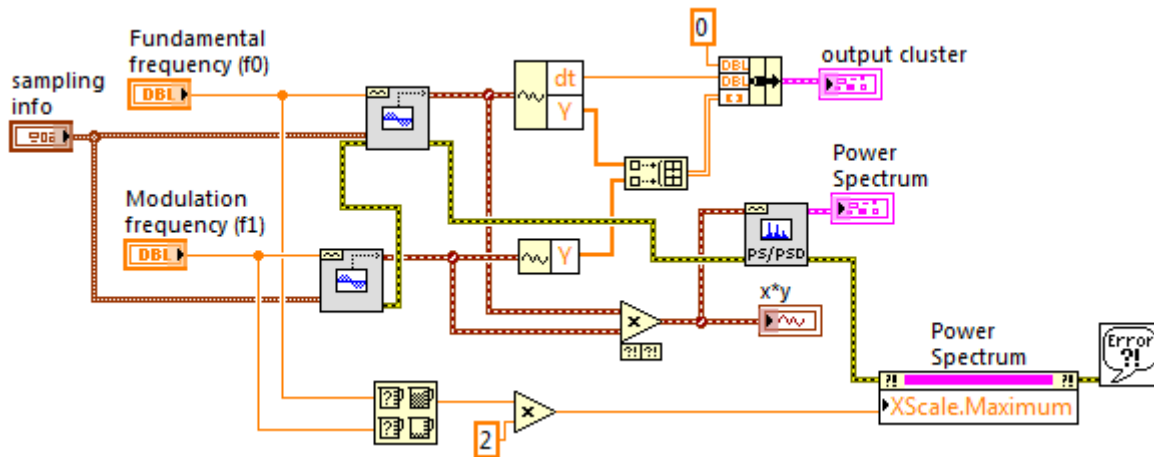


The error cluster is identified by the thick yellow-white wire. The path of the wire forces the f1 sine wave generator to execute first, followed by the f0 generator, the FFT operation, and finally the Property Node write. A Simple Error Handler VI (found in the Dialog & User Interface palette) can be optionally added to alert the user to any problems encountered in these four operations.

For the present situation, only a wire between the Power Spectrum VI error output and the Property Node error input is necessary to establish the desired sequence.


Locate spectral peaks

To increase the capability of the example VI, the frequencies of the two spectral peaks should be automatically identified. Go to the Block Diagram, right-click and select Cluster, Class, & Variant: Unbundle by Name. Wire the power spectrum bundle to its input terminal. Right-click on the icon and “Add Element”. Right-click on the top output terminal and Select Item: magnitude. Select df for the bottom terminal. Note that a 2-D array of x - and y -values is *not* available, which is why a Waveform Graph is used instead of an XY Graph. The terminal labeled “magnitude” holds a 1-D array of y -values (y_i).

Select Array: Max & Min and place it on the Block Diagram near the Unbundle operation. Wire the 1-D array from **magnitude** to the input of Min & Max. We are interested in the index of the largest y -value. The frequency corresponding to this peak is found by multiplying its index (an integer) times df . Display the result as an indicator and confirm it matches one of the peaks on the graph.

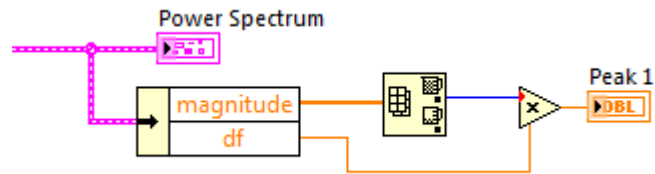

There is a second peak in the FFT spectrum that will have the same or slightly lower magnitude (this is due to sampling imprecision/noise). There are many possibilities for determining its frequency. A simple method is to replace the magnitude of the detected first peak with zero and repeating the maximum search. This is accomplished with the function Array: Replace Array Subset wired as shown below.

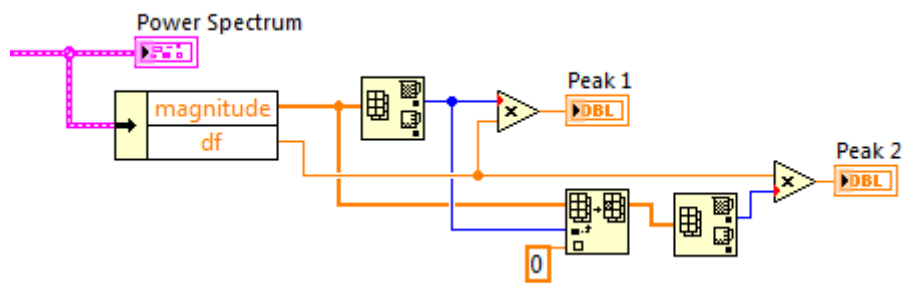

LabView has a built-in utility that can accomplish multiple spectral peak detection. Right-click on the Block Diagram and select Waveform: Analog Wfm: Measurements: Extract Multiple Tone Information VI. Wire the modulated time signal waveform to its input. Right-click on the **multiple tone information** output terminal and create an indicator on Front Panel. Run the VI and check that two frequencies are detected by scrolling through the two component cluster (select index 0 and 1 on the Front Panel).